

Analysis of Urine Biomarker for Alcohol Consumption: Screening of ETG in Urine at 9 Seconds per Sample Using LUXON-MS/MS

Pierre Picard, Jean Lacoursière and Serge Auger
Phytronix Technologies, Québec, Canada
Keywords: High-Throughput, Urine, Luxon-MS/MS

Introduction

Ethyl glucuronide is a direct metabolite of ethanol and is formed by the enzymatic conjugation of ethanol with glucuronic acid. Alcohol is normally detected for only a few hours, whereas ethyl glucuronide (ETG) can be detected in the urine for 1 to 3 days. The creatinine concentration in urine is screened to verify the sample's integrity. Normal creatinine levels ($>50 \mu\text{g/mL}$) indicate that the test sample is undiluted.

Our goal for this application note is to use an automated sample preparation method for the simultaneous quantification of creatinine (urine dilution marker) and ethyl glucuronide (ETG) in urine using a single operation in LUXON-MS/MS and perform a method comparison study according to the Trevethan¹ approach.

LUXON-MS/MS offers specificity combined with an ultra-fast analysis for an unrivaled screening method. To develop this application, we focused on performing a quick and simple sample preparation. Metabolites are analyzed **simultaneously** with **quantitative** screening results obtained in less than 9 seconds per sample. Ethyl glucuronide has been screened at a cutoff of 500 ng/mL .

Luxon Ionization Source

The Luxon Ion Source[®] (Figure 1) is the second-generation sample introduction and ionization source based on the LDTD[®] technology for mass spectrometry. Luxon Ion Source[®] uses Fiber-Coupled Laser Diode (Figure 2) to obtain unmatched thermal uniformity providing more precision, accuracy and speed. The process begins with dry samples which are rapidly evaporated using indirect heat. The thermally desorbed neutral molecules are carried into a corona discharge region. High efficiency protonation and strong resistance to ionic suppression characterize this type of ionization and is the result of the absence of solvent and mobile phase. This thermal desorption process yields high-intensity molecular ion signal in less than 1 second sample-to-sample and allows working with very small volumes.



Figure 1 - Luxon Ion Source[®]



Figure 2 - Schematic of the Luxon Ionization Source

Sample Preparation Method

Automated Sample Extraction

Urine samples were transferred into barcoded tubes, readable by the Azeo extraction system.

Each barcoded vial was scanned by the Azeo Liquid Handler and an automatic batch file was created. The Azeo extraction system (Figure 3) is used to extract the samples using the following conditions:

- 230 μL of Internal standard (ETG-d₅: 500 ng/mL and Creatinine-d₃: $4 \mu\text{g/mL}$ in dilution buffer) were added to a deep-well plate
- 20 μL of urine sample were transferred from the vials to the deep-well plate
 - Mix
- 300 μL of Dilution buffer (Methanol:Water:MTBE / 86.5:12.5:1) and 20 μL urine/internal standard mixtures were added in the deep-well plate placed on the Lumo Vortexer
 - Mix
- Spot 4 μL final dilution onto a LazWell[™]AD plate
 - Dry 4 minutes at 40°C in the Aura LazWell Dryer



Figure 3 - Automated Extraction System

LDTD[®]-MS/MS Parameters

LDTD

Model: Luxon S-960, Phytronix

Carrier gas: 6 L/min (air)

Laser pattern:

- 6-second ramp to 65% power
- Hold 2 seconds at 65% power

MS/MS

MS model: Q-Trap System[®] 5500, Sciex

Scan Time: 25 msec

Total run time: 9 seconds per sample

Ionization: APCI

Analysis Method: Negative MRM mode

Table 1 - MRM Transitions for Luxon-MS/MS

	Transition	CE
Creatinine	112 \rightarrow 68	-32
Creatinine-d ₃	115 \rightarrow 68	-32
ETG	221 \rightarrow 75	-15
ETG-d ₅	226 \rightarrow 75	-15

Results and Discussion

Initial Cut-off Test

Table 2 shows the suggested screening cut-offs currently used in the Industry.

Table 2 – Analytes and Cut-offs

Analyte	Cut-off
Ethyl glucuronide	500 ng/mL
Creatinine	50 µg/mL

Precision

Spiked samples around the decision point (50% cut-off: QC-L, cut-off: CO and 200% cut-off: QC-H) and blank solutions are used to validate the precision of the method. The peak area against the internal standard (IS) ratio was used to normalize the signal. Replicate extractions are deposited on a LazWell™ plate and dried before analysis.

The following acceptance criteria were used:

- Each concentration must not exceed 20% CV
- Mean concentration \pm 2 times the standard deviation must not overlap with other concentrations at the cut-off

For the inter-run precision experiment, each fortified sample set is analyzed in triplicate on five different days. **Table 3** shows the inter-run precision results. No overlapping at the cut-off is observed for ETG and the %CV was below 20%. Similar results are obtained for creatinine.

Table 3 - Inter-Run Precision

ETG	QC-L	Cut-off	QC-H
Conc (ng/mL)	250	500	1000
N	15	15	15
Mean (ng/mL)	247.7	529.4	967.6
SD	33.8	63.4	81.9
%CV	13.6	12.0	8.5
Mean - 2SD (ng/mL)	180.1	402.5	803.8
Mean + 2SD (ng/mL)	315.3	656.3	1131.3

For the intra-run precision experiment, each fortified sample is extracted and analyzed (8 replicates). Area ratio results are plotted using the \pm 2 STD error bars. **Figure 4** shows the intra-run results for ETG. No overlapping is observed for each concentration and the %CV was below 20%. Similar results are obtained for creatinine.

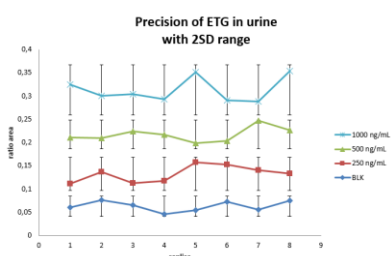


Figure 4 - Intra-Run Precision Curves for THC

Multi-Matrix Validation

Forty-five (45) real urine samples from anonymous donors were analyzed. Samples were screened to verify the presence of ETG and the endogenous level of creatinine. The drugs were analyzed with a reference method (LC-MS/MS) and the results were compared to evaluate the method sensitivity, specificity, positive predictive value (PPV), negative predictive value (NPV) and accuracy based on following calculation reported by Trevethan¹:

		LC-MS/MS	
		Yes	No
Luxon-MS/MS	Yes	TP (True positive)	FP (False positive)
	No	FN (False negative)	TN (True negative)

Where:

- Sensitivity: $(TP / (TP + FN))$
- Specificity: $(TN / (TN + FP))$
- PPV: $(TP / (TP + FP))$
- NPV: $(TN / (TN + FN))$
- Accuracy: $((TP+TN) / (TP + FN+TN+FP))$

Table 4 shows the analysis result of 45 real samples for ETG.

Table 4 – ETG Results

ETG		LC-MS/MS	
		Yes	No
Luxon-MS/MS	Yes	TP = 23	FP = 1
	No	FN = 0	TN = 21

Validation results are reported in **Table 5** for ETG.

Table 5 – Validation Results for ETG

Parameters	THC
Sensitivity (%)	100
Specificity (%)	95
PPV (%)	96
NPV (%)	100
Accuracy (%)	98

For the analysis of creatinine, all samples have a concentration greater than 50 µg/mL. Creatinine concentrations were evaluated using the isotopic analysis approach for Luxon-MS/MS and LC-MS/MS analysis. Correlation greater than 0.95 was obtained for the method comparison. **Figure 5** shows the method comparison curve.

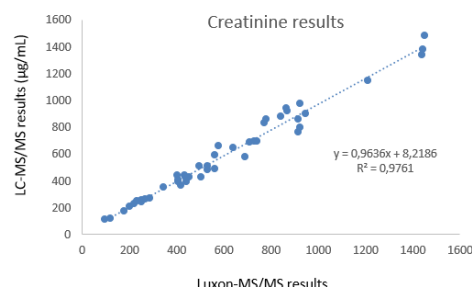


Figure 5 – Method Comparison Curve for Creatinine

Conclusion

Luxon Ion Source® combined to a Sciex Q-Trap 5500 mass spectrometer system allows ultra-fast (**9 seconds per sample**) screening of Ethyl glucuronide (ETG) and Creatinine in urine using a simple and automated sample preparation method.

Reference:

- 1) Trevethan R. Sensitivity, Specificity, and predictive values: Foundations, plabilities, and Pitfalls in research and practice. Front. Public Health (2007). 5:307

For more application notes, visit
www.phytronix.com

For research use only. Not for use in diagnostic procedures.

© 2021 Phytronix Technologies and Sciex. The trademarks mentioned herein are the property of Phytronix Technologies or Sciex or their respective owners.

Phytronix Technologies
Parc technologique du Québec métropolitain
4535, boul. Wilfrid-Hamel, Suite 120, Québec (QC)
Canada, G1P 2J7